

# Clean Circulating LV DNA Kit

Catalog Nos. CCLV-D0004, CCLV-D0096, CCLV-D0384  
 Manual revision v3.02

## Contents

Introduction and Principle .....	2
Kit Contents and Materials .....	3
Preparation of Reagents .....	5
Tube protocol for up to 1 mL Serum/Plasma .....	6
Tube protocol for up to 2 mL Serum/Plasma .....	8
Tube protocol for up to 4 mL Serum/Plasma .....	11
Plate protocol for up to 1 mL Serum/Plasma .....	14
Plate protocol for up to 2 mL Serum/Plasma .....	17
Plate protocol for up to 4 mL Serum/Plasma .....	20
Troubleshooting Guide .....	23
Ordering Information .....	25
Notes .....	26

For Research Use Only. Not for use in diagnostic procedures.  
 Information in this document is subject to change without notice.

CLEANNA DISCLAIMS ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE. TO THE FULLEST EXTENT ALLOWED BY LAW, IN NO EVENT SHALL CLEANNA BE LIABLE, WHETHER CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING, BUT NOT LIMITED TO THE USE THEREOF, WHETHER OR NOT FORSEEABLE AND WHETHER OR NOT CLEANNA IS ADVISED OF THE POSSIBILITY OF SUCH DAMAGES.



## Introduction and Principle

The Clean Circulating LV DNA Kit allows for the isolation of cell free DNA from sample volumes up to 4 mL plasma or serum. The entire procedure allows the kit to be used both manually as well as automated.

Its unique buffer system, eliminates the needs for vacuum steps as well as funnels throughout the procedure. As a result, the Clean Circulating LV DNA Kit provides hands-free procedure for automated protocols via a simple 4 step process (lyse, bind, wash and elute).

Our CleanNA Particles CCLV offer a high binding capacity and combined with our propriety buffer system target smaller DNA fragments (120-400 bp). This combination minimizes the risk of genomic DNA contamination.

The high binding capacity of the CleanNA particles CCLV decrease the amount of particles required during binding steps thereby reducing the elution volume. This enables isolated cell free DNA from 4 mL plasma to be eluted in as little as 50  $\mu$ L.

Using a special formulated lysis buffer, the circulating DNA is released from proteins or vesicles bound to the DNA while DNases and RNases are inactivated. DNA is isolated from the lysates in one step by binding to CleanNA Particles' surfaces. The CleanNA magnetic particles are separated from the lysates by using a magnetic separation device. Following a few rapid wash steps to remove trace contaminants, the purified DNA is eluted from the CleanNA particles for downstream applications using an Elution Buffer.

The isolated Cell Free DNA is directly suitable for use in downstream applications such as Next Generation Sequencing (NGS) and qPCR.

## Kit Contents and Materials

### Kit Contents:

Product	CCLV-D0004	CCLV-D0096	CCLV-D0384	Storage
Preps	4	96	384	
CleanNA Particles CCLV	200 µL	4 mL	13 mL	2-8°C
CCLV Lysis	1.5 mL	30 mL	120 mL	20-25°C
CCLV Binding	20 mL	450 mL	2 x 850 mL	20-25°C
CCLV Wash 1	10 mL	2 x 120 mL	900 mL	20-25°C
CCLV Wash 2	2.5 mL	3 x 20 mL	200 mL	20-25°C
Elution Buffer	10 mL	250 mL	1000 mL	20-25°C
Proteinase K Solution	300 µL	7 mL	28 mL	20-25°C (for storage > 12 months, store at 2-8°C)

### Materials to be supplied by user:

#### For isolation in single tubes:

#### Materials and Reagents to be supplied by User for the Tube Protocol for up to 1 mL of sample input:

- 100% ethanol
- Magnetic separation device for 1.5/2.0 mL tubes
- Vortexer
- Shaker or Rocker
- Incubator capable of 60°C
- 1.5 mL micro centrifuge tube(s)
- 15 mL centrifuge tube(s)

#### Materials and Reagents to be supplied by User for the Tube Protocol for up to 2 mL and up to 4 mL of sample input:

- 100% ethanol
- Magnetic separation device for 15 mL centrifuge tubes
- Magnetic separation device for 1.5 / 2.0 mL tubes
- Vortexer
- Shaker or Rocker
- Incubator capable of 60°C
- 1.5 mL micro centrifuge tube(s)
- 15 mL centrifuge tube(s), compatible with magnetic separation device used

**For isolation using 96-well plate format:**

**Materials and Reagents to be supplied by User for the Plate Protocol for up to 1 mL of sample input:**

- 100% ethanol
- Magnetic separation device for 96-well plates from CleanNA (Cat# CMAG-RN50)
- Vortexer
- Shaker or Rocker
- Incubator capable of 60°C
- 24-well deep-well plate(s), 10 mL (Cat# Whatman 7701-5102)
- 96-well deep-well plate(s)

**Materials and Reagents to be supplied by User for the Plate Protocol for up to 2 mL and up to 4 mL of sample input:**

- 100% ethanol
- Magnetic separation device for 96-well plates from CleanNA (Cat# CMAG-RN50)
- Vortexer
- Shaker or Rocker
- Incubator capable of 60°C
- 24-well deep-well plate(s), 10 mL (Cat# Whatman 7701-5102)
- 96-well deep-well plate(s)

## Preparation of Reagents

Dilute CCLV Wash 2 with 100% ethanol as follows and store at room temperature.

Kit	100% Ethanol to be Added
CCLV-D0004	10 mL
CCLV-D0096	80 mL
CCLV-D0384	800 mL

## Clean Circulating LV Kit

### Tube protocol for up to 1 mL Serum/Plasma

#### Before Starting:

- Set Incubator to 60°C
- Prepare CCLV Wash 2 according to the instructions in the Preparing Reagents section on page 5.
- Shake or Vortex the CleanNA Particles CCLV to fully resuspend the particles before use.

#### Protocol:

1. Add up to 1 mL plasma/serum samples to a 15 mL centrifuge tube (not provided).
2. Bring the sample volume up to 1 mL with Elution Buffer (provided with this kit) if the sample volume is less than 1 mL.
3. Add 15 µL Proteinase K Solution.
4. Add 67 µL CCLV Lysis.
5. Vortex at maximum speed or pipet up and down to mix thoroughly.
6. Incubate at 60°C for 20 minutes. Mix by inverting or shaking every 10 minutes.
7. Incubate at room temperature for 10 minutes.





**Note:** This incubation step is crucial to let the sample temperature drop and obtain most efficient DNA binding to the CleanNA Particles CCLV.

8. Add 1 mL CCLV Binding. Vortex at maximum speed for 30 seconds or pipet up and down to mix thoroughly.
9. Add 10 µL CleanNA Particles CCLV. Invert the sample 10 times or pipet up and down to mix.
10. Incubate for 10 minutes at room temperature with continuous mixing. The samples must be mixed throughout the 10 minute incubation period by shaking or rocking.



**Note:** Do not vortex at high speeds as this will cause foaming resulting in a reduced yield. The speed of mixing should be set to continuously keep the CleanNA Particles CCLV resuspended in solution.

11. Transfer 1 mL lysate to a 1.5 mL micro centrifuge tube (not provided).
12. Place the tube on a magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
13. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
14. Transfer the remaining lysate from step 11 to the 1.5 mL micro centrifuge tube used in the previous steps.
15. Place the tube on a magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
16. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
17. Remove the tube containing the CleanNA Particles CCLV from the magnetic separation device.

18. Add 500 µL CCLV Wash 1.
19. Resuspend the CleanNA Particles CCLV by vortexing for 2 minutes or pipetting up and down 20 times.
  -  **Note:** To obtain good purity, complete resuspension of the CleanNA Particles CCLV is critical.
20. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
21. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
22. Repeat steps 17-21 for a second “CCLV Wash 1” wash step
23. Remove the tube containing the CleanNA Particles CCLV from the magnetic separation device.
24. Add 500 µL CCLV Wash 2.
  -  **Note:** CCLV Wash 2 must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.
25. Resuspend the CleanNA Particles CCLV by vortexing for 2 minutes or pipetting up and down 20 times.
26. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
27. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
28. Repeat Steps 23-27 for a second “CCLV Wash 2” wash step.
29. Remove the tube from the magnetic separation device for approximately 30 seconds.
30. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV.
31. Aspirate and discard the residual CCLV Wash 2.
32. Leave the tube on the magnetic separation device for 25 minutes to dry the CleanNA Particles CCLV.
33. Remove the tube containing the CleanNA Particles CCLV from the magnetic separation device.
34. Add 30-60 µL Elution Buffer. Resuspend the CleanNA Particles CCLV by vortexing or pipetting up and down 20 times.
35. Incubate at room temperature for 5 minutes, while constantly vortexing.
36. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
37. Transfer the cleared supernatant containing purified DNA to a clean 1.5 mL micro centrifuge tube (not provided).
38. Store DNA at -20°C.

## Clean Circulating LV Kit Tube protocol for up to 2 mL Serum/Plasma

### Before Starting:

- Set Incubator to 60°C
- Prepare CCLV Wash 2 according to the instructions in the Preparing Reagents section on page 5.
- Shake or Vortex the CleanNA Particles CCLV to fully resuspend the particles before use.

### Protocol:

1. Add up to 2 mL plasma/serum samples to a 15 mL centrifuge tube (not provided). Choose the correct plastic ware depending on the magnetic separation device being used.
2. Bring the sample volume up to 2 mL with Elution Buffer (provided with this kit) if the sample volume is less than 2 mL.
3. Add 30 µL Proteinase K Solution.
4. Add 135 µL CCLV Lysis.
5. Vortex at maximum speed or pipet up and down to mix thoroughly.
6. Incubate at 60°C for 25 minutes. Mix by inverting or shaking every 10 minutes.
7. Incubate at room temperature for 10 minutes.



**Note:** This incubation step is crucial to let the sample temperature drop and obtain most efficient DNA binding to the CleanNA Particles CCLV.

8. Add 2 mL CCLV Binding. Vortex at maximum speed for 30 seconds or pipet up and down to mix thoroughly.
9. Add 20 µL CleanNA Particles CCLV. Invert the sample 10 times or pipet up and down to mix.
10. Incubate for 10 minutes at room temperature with continuous mixing. The samples must be mixed throughout the 10 minute incubation period by shaking or rocking.



**Note:** Do not vortex at high speeds as this will cause foaming resulting in a reduced yield. The speed of mixing should be set to continuously keep the CleanNA Particles CCLV resuspended in solution.

11. Place the tube on a magnetic separation device, compatible with the 15 mL tube, to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
12. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
13. Remove the tube/plate containing the CleanNA Particles CCLV from the magnetic separation device.
14. Add 1 mL CCLV Wash 1.
15. Resuspend the CleanNA Particles CCLV by vortexing for 2 minutes or pipetting up and down 20 times.



**Note:** To obtain good purity, complete resuspension of the CleanNA Particles CCLV is critical.



16. Transfer the resuspended CleanNA Particles CCLV to a new 1.5 mL microcentrifuge tube (not provided).



**Note:** Use a magnetic separation device designed for 1.5/2.0 mL tubes for the remaining procedure.

17. Place the 1,5 mL tube on the magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
18. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
19. Remove the tube containing the CleanNA Particles CCLV from the magnetic separation device.
20. Add 1 mL CCLV Wash 1.
21. Resuspend the CleanNA Particles CCLV by vortexing for 2 minutes or pipetting up and down 20 times.



**Note:** To obtain good purity, complete resuspension of the CleanNA Particles CCLV is critical.

22. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
23. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
24. Remove the tube containing the CleanNA Particles CCLV from the magnetic separation device.
25. Add 1 mL CCLV Wash 2.



**Note:** CCLV Wash 2 must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

26. Resuspend the CleanNA Particles CCLV by vortexing for 2 minutes or pipetting up and down 20 times.
27. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
28. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
29. Repeat Steps 24-28 for a second CCLV Wash 2 wash step.
30. Remove the tube from the magnetic separation device for approximately 30 seconds.
31. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV.
32. Aspirate and discard the residual CCLV Wash 2.
33. Leave the tube on the magnetic separation device for 25 minutes to dry the CleanNA Particles CCLV.
34. Remove the tube containing the CleanNA Particles CCLV from the magnetic separation device.
35. Add 50-100  $\mu$ L Elution Buffer. Resuspend the CleanNA Particles CCLV by vortexing or pipetting up and down 20 times.

36. Incubate at room temperature for 5 minutes, while constantly vortexing.
37. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
38. Transfer the cleared supernatant containing purified DNA to a clean 1.5 mL micro centrifuge tube (not provided).
39. Store DNA at -20°C.

## Clean Circulating LV Kit

### Tube protocol for up to 4 mL Serum/Plasma

#### Before Starting:

- Set Incubator to 60°C
- Prepare CCLV Wash 2 according to the instructions in the Preparing Reagents section on page 5.
- Shake or Vortex the CleanNA Particles CCLV to fully resuspend the particles before use

#### Protocol:

1. Add up to 4 mL plasma/serum samples to a 15 mL centrifuge tube (not provided). Choose the correct plastic ware depending on the magnetic separation device being used.
2. Bring the volume up to 4 mL with Elution Buffer (provided with this kit) if the sample volume is less than 4 mL.
3. Add 60 µL Proteinase K Solution.
4. Add 270 µL CCLV Lysis.
5. Vortex at maximum speed or pipet up and down to mix thoroughly.
6. Incubate at 60°C for 30 minutes. Mix by inverting or shaking every 10 minutes.
7. Incubate at room temperature for 10 minutes.



**Note:** This incubation step is crucial to let the sample temperature drop and obtain most efficient DNA binding to the CleanNA Particles CCLV.

8. Add 4 mL CCLV Binding. Vortex at maximum speed for 30 seconds or pipet up and down to mix thoroughly.
9. Add 30 µL CleanNA Particles CCLV. Invert the sample 10 times or pipet up and down to mix.
10. Incubate for 10 minutes at room temperature with continuous mixing. The samples must be mixed throughout the 10 minute incubation period by shaking or rocking.



**Note:** Do not vortex at high speeds as this will cause foaming resulting in a reduced yield. The speed of mixing should be set to continuously keep the CleanNA Particles CCLV resuspended in solution.

11. Place the tube on a magnetic separation device, compatible with the 15 mL tube, to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
12. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
13. Remove the tube containing the CleanNA Particles CCLV from the magnetic separation device.
14. Add 1 mL CCLV Wash 1.
15. Resuspend the CleanNA Particles CCLV by vortexing for 5 minutes or pipetting up and down 20 times.



**Note:** To obtain good purity, complete resuspension of the CleanNA Particles CCLV is critical.

16. Transfer the resuspended CleanNA Particles CCLV to a new 1.5 mL micro centrifuge tube (not provided).



**Note:** Use a magnetic separation device designed for 1.5/2.0 mL tubes for the remaining procedure.

17. Place the 1.5 mL tube on the magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.

18. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.

19. Remove the tube containing the CleanNA Particles CCLV from the magnetic separation device.

20. Add 1 mL CCLV Wash 1.

21. Resuspend the CleanNA Particles CCLV by vortexing for 5 minute or pipetting up and down 20 times.



**Note:** To obtain good purity, complete resuspension of the CleanNA Particles CCLV is critical.

22. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.

23. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.

24. Remove the tube containing the CleanNA Particles CCLV from the magnetic separation device.

25. Add 1 mL CCLV Wash 2.



**Note:** CCLV Wash 2 must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

26. Resuspend the CleanNA Particles CCLV by vortexing for 5 minutes or pipetting up and down 20 times.

27. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.

28. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.

29. Repeat Steps 24-28 for a second CCLV Wash 2 wash step.

30. Remove the tube from the magnetic separation device for approximately 30 seconds.

31. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV.

32. Aspirate and discard the residual CCLV Wash 2.

33. Leave the tube on the magnetic separation device for 25 minutes to dry the CleanNA Particles CCLV.

34. Remove the tube containing the CleanNA Particles CCLV from the magnetic separation device.

35. Add 50-100  $\mu$ L Elution Buffer. Resuspend the CleanNA Particles CCLV by vortexing or pipetting up and down 20 times.

36. Incubate at room temperature for 5 minutes, while constantly vortexing.
37. Place the tube on the magnetic separation device to magnetize the CleanNA Particles CCLV. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
38. Transfer the cleared supernatant containing purified DNA to a clean microplate or 1.5 mL micro centrifuge tube (not provided).
39. Store DNA at -20°C.

## Clean Circulating LV Kit

### Plate protocol for up to 1 mL Serum/Plasma

#### Before Starting:

- Set Incubator to 60°C
- Prepare CCLV Wash 2 according to the instructions in the Preparing Reagents section on page 5.
- Shake or Vortex the CleanNA Particles CCLV to fully resuspend the particles before use.

#### Protocol:

1. Add up to 1 mL plasma/serum samples to a 24-well Deep Well Plate (not provided).
2. Bring the sample volume up to 1 mL with Elution Buffer (provided with this kit) if the sample volume is less than 1 mL.
3. Add 15 µL Proteinase K Solution.
4. Add 67 µL CCLV Lysis, optionally seal the plate.
5. Vortex at maximum speed or pipet up and down to mix thoroughly.
6. Incubate at 60°C for 20 minutes. Mix by inverting or shaking every 10 minutes.
7. Incubate at room temperature for 10 minutes.



**Note:** This incubation step is crucial to let the sample temperature drop and obtain most efficient DNA binding to the CleanNA Particles CCLV.

8. Add 1 mL CCLV Binding. Vortex at maximum speed for 30 seconds or pipet up and down to mix thoroughly.
9. Add 10 µL CleanNA Particles CCLV. Invert the sample 10 times or pipet up and down to mix.
10. Incubate for 10 minutes at room temperature with continuous mixing. The samples must be mixed throughout the 10 minute incubation period by shaking or rocking.



**Note:** Do not vortex at high speeds as this will cause foaming resulting in a reduced yield. The speed of mixing should be set to continuously keep the CleanNA Particles CCLV resuspended in solution.

11. Place the 24-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV. The Particles from each well will be collected by for magnets at the bottom.
12. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
13. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
14. Remove the 24-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
15. Add 500 µL CCLV Wash 1.
16. Resuspend the CleanNA Particles CCLV by vortexing for 2 minutes or pipetting up and down 20 times.



**Note:** To obtain good purity, complete resuspension of the CleanNA Particles CCLV is critical.

17. Transfer the resuspended CleanNA Particles CCLV to a new 96-well Deep Well Plate (not provided).



**Note:** Continue to work in 96-well format for the remaining procedure.

18. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
19. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
20. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
21. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
22. Add 500  $\mu$ L CCLV Wash 1.

23. Resuspend the CleanNA Particles CCLV by vortexing for 2 minutes or pipetting up and down 20 times.



**Note:** To obtain good purity, complete resuspension of the CleanNA Particles CCLV is critical.

24. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
25. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.  
  
Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
26. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
27. Add 500  $\mu$ L CCLV Wash 2.



**Note:** CCLV Wash 2 must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

28. Resuspend the CleanNA Particles CCLV by vortexing for 2 minutes or pipetting up and down 20 times.
29. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
30. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
31. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
32. Repeat Steps 26-31 for a second “CCLV Wash 2” wash step.
33. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
34. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
35. Aspirate and discard the residual CCLV Wash 2.

36. Leave the tube on the magnetic separation device for 25 minutes to dry the CleanNA Particles CCLV.
37. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
38. Add 30-60  $\mu$ L Elution Buffer. Resuspend the CleanNA Particles CCLV by vortexing or pipetting up and down 20 times.
39. Incubate at room temperature for 5 minutes, while constantly vortexing.
40. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
41. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
42. Transfer the cleared supernatant containing purified DNA to a clean 96-well plate of to clean individual tubes (not provided).
43. Store DNA at  $-20^{\circ}\text{C}$ .



## Clean Circulating LV Kit

### Plate protocol for up to 2 mL Serum/Plasma

#### Before Starting:

- Set Incubator to 60°C
- Prepare CCLV Wash 2 according to the instructions in the Preparing Reagents section on page 5.
- Shake or Vortex the CleanNA Particles CCLV to fully resuspend the particles before use.

#### Protocol:

1. Add up to 2 mL plasma/serum samples to a 24-well Deep Well Plate (not provided).
2. Bring the sample volume up to 2 mL with Elution Buffer (provided with this kit) if the sample volume is less than 2 mL.
3. Add 30 µL Proteinase K Solution.
4. Add 135 µL CCLV Lysis, optionally seal the plate.
5. Vortex at maximum speed or pipet up and down to mix thoroughly.
6. Incubate at 60°C for 25 minutes. Mix by inverting or shaking every 10 minutes.
7. Incubate at room temperature for 10 minutes.



**Note:** This incubation step is crucial to let the sample temperature drop and obtain most efficient DNA binding to the CleanNA Particles CCLV.

8. Add 2 mL CCLV Binding. Vortex at maximum speed for 30 seconds or pipet up and down to mix thoroughly.
9. Add 20 µL CleanNA Particles CCLV. Invert the sample 10 times or pipet up and down to mix.
10. Incubate for 10 minutes at room temperature with continuous mixing. The samples must be mixed throughout the 10 minute incubation period by shaking or rocking.



**Note:** Do not vortex at high speeds as this will cause foaming resulting in a reduced yield. The speed of mixing should be set to continuously keep the CleanNA Particles CCLV resuspended in solution.

11. Place the 24-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV. The Particles from each well will be collected by for magnets at the bottom.
12. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
13. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
14. Remove the 24-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
15. Add 1 mL CCLV Wash 1.
16. Resuspend the CleanNA Particles CCLV by vortexing for 2 minutes or pipetting up and down 20 times.



**Note:** To obtain good purity, complete resuspension of the CleanNA Particles CCLV is critical.

17. Transfer the resuspended CleanNA Particles CCLV to a new 96-well Deep Well Plate (not provided).



**Note:** Continue to work in 96-well format for the remaining procedure.

18. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
19. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
20. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
21. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
22. Add 1 mL CCLV Wash 1.

23. Resuspend the CleanNA Particles CCLV by vortexing for 2 minutes or pipetting up and down 20 times.



**Note:** To obtain good purity, complete resuspension of the CleanNA Particles CCLV is critical.

24. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
25. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.  
  
Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
26. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
27. Add 1 mL CCLV Wash 2.



**Note:** CCLV Wash 2 must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

28. Resuspend the CleanNA Particles CCLV by vortexing for 2 minutes or pipetting up and down 20 times.
29. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
30. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
31. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
32. Repeat Steps 26-31 for a second “CCLV Wash 2” wash step.
33. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
34. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
35. Aspirate and discard the residual CCLV Wash 2.

36. Leave the tube on the magnetic separation device for 25 minutes to dry the CleanNA Particles CCLV.
37. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
38. Add 50-100  $\mu$ L Elution Buffer. Resuspend the CleanNA Particles CCLV by vortexing or pipetting up and down 20 times.
39. Incubate at room temperature for 5 minutes, while constantly vortexing.
40. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
41. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
42. Transfer the cleared supernatant containing purified DNA to a clean 96-well plate of to clean individual tubes (not provided).
43. Store DNA at  $-20^{\circ}\text{C}$ .

## Clean Circulating LV Kit

### Plate protocol for up to 4 mL Serum/Plasma

#### Before Starting:

- Set Incubator to 60°C
- Prepare CCLV Wash 2 according to the instructions in the Preparing Reagents section on page 5.
- Shake or Vortex the CleanNA Particles CCLV to fully resuspend the particles before use.

#### Protocol:

1. Add up to 4 mL plasma/serum samples to a 24-well Deep Well Plate (not provided).
2. Bring the sample volume up to 4 mL with Elution Buffer (provided with this kit) if the sample volume is less than 4 mL.
3. Add 60 µL Proteinase K Solution.
4. Add 270 µL CCLV Lysis, optionally seal the plate.
5. Vortex at maximum speed or pipet up and down to mix thoroughly.
6. Incubate at 60°C for 30 minutes. Mix by inverting or shaking every 10 minutes.
7. Incubate at room temperature for 10 minutes.



**Note:** This incubation step is crucial to let the sample temperature drop and obtain most efficient DNA binding to the CleanNA Particles CCLV.

8. Add 4 mL CCLV Binding. Vortex at maximum speed for 30 seconds or pipet up and down to mix thoroughly.
9. Add 30 µL CleanNA Particles CCLV. Invert the sample 10 times or pipet up and down to mix.
10. Incubate for 10 minutes at room temperature with continuous mixing. The samples must be mixed throughout the 10 minute incubation period by shaking or rocking.



**Note:** Do not vortex at high speeds as this will cause foaming resulting in a reduced yield. The speed of mixing should be set to continuously keep the CleanNA Particles CCLV resuspended in solution.

11. Place the 24-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV. The Particles from each well will be collected by for magnets at the bottom.
12. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
13. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
14. Remove the 24-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
15. Add 1 mL CCLV Wash 1.
16. Resuspend the CleanNA Particles CCLV by vortexing for 5 minutes or pipetting up and down 20 times.



**Note:** To obtain good purity, complete resuspension of the CleanNA Particles CCLV is critical.

17. Transfer the resuspended CleanNA Particles CCLV to a new 96-well Deep Well Plate (not provided).



**Note:** Continue to work in 96-well format for the remaining procedure.

18. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
19. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
20. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
21. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
22. Add 1 mL CCLV Wash 1.

23. Resuspend the CleanNA Particles CCLV by vortexing for 5 minutes or pipetting up and down 20 times.



**Note:** To obtain good purity, complete resuspension of the CleanNA Particles CCLV is critical.

24. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
25. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.  
  
Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
26. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
27. Add 1 mL CCLV Wash 2.



**Note:** CCLV Wash 2 must be diluted with 100% ethanol prior to use. Please see Page 5 for instructions.

28. Resuspend the CleanNA Particles CCLV by vortexing for 5 minutes or pipetting up and down 20 times.
29. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
30. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
31. Aspirate and discard the cleared supernatant. Do not disturb the CleanNA Particles CCLV.
32. Repeat Steps 26-31 for a second “CCLV Wash 2” wash step.
33. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
34. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
35. Aspirate and discard the residual CCLV Wash 2.

36. Leave the tube on the magnetic separation device for 25 minutes to dry the CleanNA Particles CCLV.
37. Remove the 96-well plate containing the CleanNA Particles CCLV from the magnetic separation device.
38. Add 50-100  $\mu$ L Elution Buffer. Resuspend the CleanNA Particles CCLV by vortexing or pipetting up and down 20 times.
39. Incubate at room temperature for 5 minutes, while constantly vortexing.
40. Place the 96-well Deep Well Plate on the 96 ring magnet plate to magnetize the CleanNA Particles CCLV.
41. Incubate at room temperature until the CleanNA Particles CCLV are completely cleared from solution.
42. Transfer the cleared supernatant containing purified DNA to a clean 96-well plate of to clean individual tubes (not provided).
43. Store DNA at  $-20^{\circ}\text{C}$ .

## Troubleshooting Guide

Please use this guide to troubleshoot any problems that may arise. For further assistance, please contact your local distributor.

### Possible Problems and Suggestions

Problem	Cause	Solution
Low DNA yield	Incomplete resuspension of CleanNA Particles CCLV	Resuspend the CleanNA Particles CCLV by vortexing vigorously before use
	Inefficient binding of the DNA to the CleanNA Particles CCLV	Ensure to let the sample cool at room temperature for 10 minutes prior to addition of the CCLV Binding.
	Inefficient binding of the DNA to the CleanNA Particles CCLV	Ensure to mix each sample continuously throughout the binding incubation
	Loss of CleanNA Particles CCLV during operation	Avoid disturbing the CleanNA Particles CCLV during aspiration
	DNA remains bound to CleanNA Particles CCLV	Increase elution volume and incubate at room temperature for 15 minutes; Pipet up and down 50 to 100 times.
	DNA washed off	Dilute CCLV Wash 2 by adding appropriate volume of 100% ethanol prior to use (see page 5 for instructions).
	Ethanol carryover	Dry the CleanNA Particles CCLV at room temperature for 25 minutes before elution.
CleanNA particles CCLV do not completely clear from solution	Too short of magnetizing time	Increase collection time on the magnetic separation device
High Molecular Weight Co-Purification	Two CCLV Wash 1 Steps must be performed	Perform two CCLV Wash 1 steps as instructed in the manual. Increase the volume of wash buffer if necessary.
Problems in downstream applications	Salt carryover	CCLV Wash 2 must be at room temperature.

Problem	Cause	Solution
Abnormal BioAnalyzer data	BioAnalyzer shows multiple sharp peaks during analysis	Ensure to remove all traces of the cleared supernatant after each wash step
		Ensure to incubate the tube/plate for 25 minutes to dry the CleanNA Particles CCLV
	BioAnalyzer shows base line climbing towards the end	Check the BioAnalyzer chip for air bubbles. Load samples onto a new freshly prepared chip
	BioAnalyzer shows high blob at the beginning of the trace	Ensure the purified sample does not contain traces of CleanNA particles CCLV



## Ordering Information

Contact your local distributor to order.

Product	Part Number
Clean Circulating LV DNA Kit (4 Preps)	CCLV-D0004
Clean Circulating LV DNA Kit (96 Preps)	CCLV-D0096
Clean Circulating LV DNA Kit (384 Preps)	CCLV-D0384

Product	Part Number
Clean Magnet Plate 96-Well	CMAG-96-RN50

**Notes**

**Notes**

CLEAN  NA